DNA EXTRACTION FROM FRENCH SOILS EVALUATION OF PRECELLYS® 24-DUAL 7ML VIAL

Genosol – UMR Microbiology of the soil and environment, Inra Dijon, France

/ CONTEXT

The Genosol platform is a biological resource centre unique in France and Europe, devoted to the conservation and analysis of the genetic resources of soil microbial communities. The ambition of the GenoSol platform is to centralize the microbial genetic resources of soils and make them available to the scientific community together with the most up-to-date technologies for their characterization. In this way, the **Precellys® 24-Dual** was evaluated in comparison with the reference method.

/ MATERIALS

- **Precellys® 24-Dual** with Precellys® kit 7mL vial with 0.1 mm glass and 5.0 mm ceramics beads (P000944-LYSK0-A.0, P000929-LYSK0-A.0 and P000930-LYSK0-A.0).
- Standard method: Mikro-Dismembrator S (Sartorius).
- Samples: Three clayey, silty-clayey, sandy soils (Arceau, Champdôtte, Epoisses).

/ PROTOCOL


- **Precellys® 24-Dual** protocol : 1g soil + 4mL Buffer / 7mL vial, 5000 rpm 3x15s. [alternatively, on **Precellys Evolution**, the following program can be used: 7200rpm 3x15s]
- Electrophoresis gel
- DNA extraction yield, qPCR.

/ CONCLUSION


The **Precellys® 24-Dual** is successfully validated for extraction of total DNA with higher yield of extraction. The DNA quality is optimal to study the soil microbial communities. Easy handling, reproducibility and high throughput are appreciated.

/ RESULTS

With **Precellys® 24-Dual**, the concentration of DNA extracted is around 2.5 higher than the standard method. The quality of DNA, evaluated by inhibition and restrictive enzyme assays was equivalent with both equipments. **Precellys® 24-Dual** allows a better cadence: 6 samples could be prepared in 45s versus 120s with standard method (1 sample by run).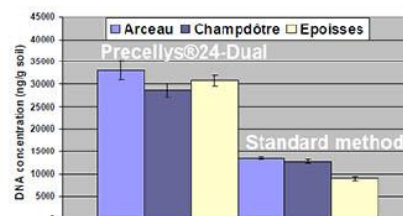

DNA concentration (ng/g soil) for 3 French soils.

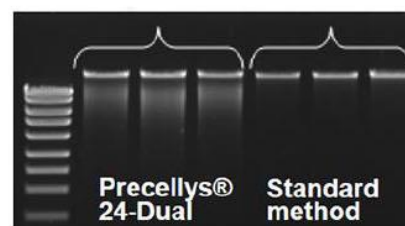Gel electrophoresis analysis, Epoisses soil (triplicate).

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